

### Safety Instructions



Wear safety goggles at all times.

Acids and bases are corrosive. Do not get any on skin or near eyes.

In case of a spill, rinse with large amounts of water.

### pH Paper Test Procedure

1. Place your pH paper on the watch glass in front of the solution.
2. Use the glass stirring rod to place one drop of solution on the end of the pH paper.
3. Record your observations in the table.
4. Discard your pH paper in the waste container. Rinse the watch glass.

### pH Meter Test Procedure

1. Rinse the probe with distilled water.
2. Place the probe in the solution to be tested.
3. Record the pH in your data table.
4. Rinse the probe with distilled water before moving to the next solution.
5. When finished, rinse the probe with distilled water and place it in the storage beaker of distilled water.